DOSING  
SYSTEM

CO<sub>2</sub>RA<sup>®</sup>

# DOSING SYSTEM

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*Every process of the chemical, pharmaceutical, food and homeopathic industry uses powders, granules, tablets and capsules and requires dosing systems for the intermediate packaging of the product.*

*CO.RA. offers customized solutions suited to meet the needs and solve the specific problems due to the type of product or process where a dosing system is necessary. The DOSING SYSTEM allows you to discharge and dose products such as powders, granules, tablets and capsules in bags, containers or drums which must be stored temporarily before being sent to the end-user and where a maximum weight accuracy could be required.*

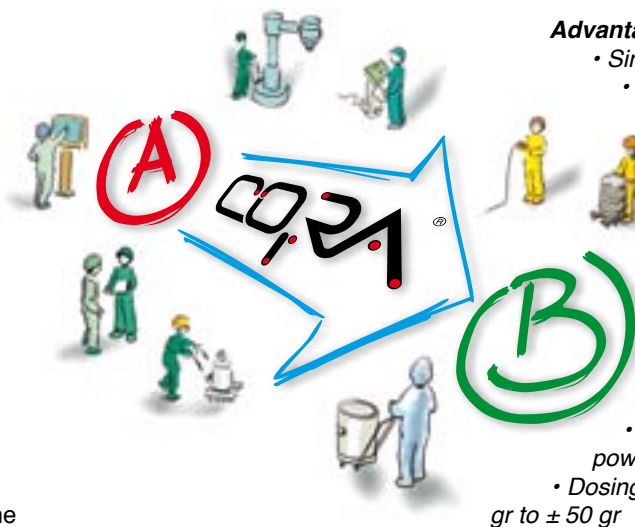
*The DOSING SYSTEM comes in four different versions:*

- 1 Inflatable ring dosing system
  - 2 Continuous liner dosing system
  - 3 Dosing system with "Y"
  - 4 Dosing system for tablets and capsules
- and these four solutions can be produced according to the end-user's needs.*

*The thirty years of experience in dealing with the complexity stringent standards governing the above mentioned sectors, in particular the pharmaceuticals, and our choice to deliver products in AISI 316/304 has allowed us to introduce dedicated machines for the discharging, dosing and handling of powders to meet the needs of the chemical, pharmaceutical, homeopathic and food industries.*

*The stainless steel parts not in contact with the product have a satin finish (Ra 0,8) and the contact parts are mirror polished (Ra 0.2). On request and based on the area in which it will be installed, the dosing system can be entirely mirror polished.*

*CO.RA. dosing systems are easy to integrate within existing production processes, but may also be part of a "turnkey" project to meet the customer's specific requirements.*

## Advantages:

- Simple and compact design
- Designed and manufactured in accordance to the end-user's specific requirements
- Easy to disassemble, inspect and clean
- Materials: AISI 304 and 316L with silicone seals (certificates on request)
- Manufactured according to GMP norms (Good Manufacturing Practice)
- Quick and easy to maintenance
- Adaptable for different types of powders, capsules and tablets
- Dosing valves with an accuracy from  $\pm 500$  gr to  $\pm 50$  gr

## Applications:

- Where it is necessary to dose into drums or bags
- Weighing and filling systems for bags, containers and drums
- Intermediate stocking

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### 1. INFLATABLE RING DOSING SYSTEM

*The inflatable ring dosing system is used in processes where the end-user doses the product inside of bags transported in containers, where a dosing accuracy between  $\pm 500$ g and  $\pm 50$ g is required. This system not only ensures a reduced contact with the product, but it is also easy to use and clean. The main feature of this system is the inflatable ring, which can be supplied in two versions:*

**Inflatable:** when in the resting position the ring is deflated, but when the work cycle begins the ring inflates and blocks the sack against the frame, preventing the bag from being pulled away by the weight of the product being discharged.

**Deflatable:** when in the resting position the ring is inflated, but when the work cycle begins, it deflates creating the necessary space for the operator to insert the bag. Once the bag has been inserted the ring is inflated again. As for the "Inflatable" version, the ring blocks the sack against the frame, preventing the bag from being pulled out by the weight of the product being discharged.

The system is composed of:

- Interface adaptor for the machine to be served
- Discharge pipeline
- Intercepting valve which can be of two different types:
  - DOUBLE VALVE and SANITARY ROTARY VALVE for a dosing accuracy between  $\pm 100$  gr and  $50$  gr
  - ROTARY VALVE for bulk dosing where required accuracy is between  $\pm 500$  gr and  $100$  gr.
- Inflatable ring
- Supporting structure which can be fixed and/or on wheels
- Weighing system
- Control panel

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## 2. CONTINUOUS LINER DOSING SYSTEMS

*The continuous liner dosing system is used in processes where the end-user doses products in drums, in absence of a bag. This type of system is used in processes where the end-user doses the product inside of bags transported in containers, where the dosing accuracy required is between  $\pm 500$ g and  $\pm 50$ g.*

*The operator puts the container on the scale, closes the bag, extends the liner to the requested height and then begins to discharge. Once it is finished the operator closes the bag again. The methods of closing the bag are two: with a disposable clip or with a sealer and in both cases this type of system allows the operator to close the bag avoiding the leakage of the product into the work environment and ensuring the reduced contact of the operator with the product.*

*This system is composed of:*

- Interface adaptor for the machine to be served
- Discharge pipeline
- Intercepting valve which can be of two different types:
  - DOUBLE VALVE and SANITARY ROTARY VALVE for a dosing accuracy between  $\pm 100$  gr and 50 gr
  - ROTARY VALVE for bulk dosing where required accuracy is between  $\pm 500$  gr and 100 gr.
- Liner holding structure
- Supporting structure which can be fixed and/or on wheels
- Weighing system
- Control panel

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### 3. DOSING SYSTEM WITH "Y"

*The dosing system with "Y" is a system used in processes where the product must be discharged into two drums, or in a drum and a bin at the same time. The "Y" pipeline is installed directly in its working position and depending on the type of product and the type of dosing required, a ROTARY VALVE or VR8V can be installed on the inlet.*

*The ROTARY VALVE allows the discharge of large quantities of product by means of a reciprocating 180° movement, the ROTARY VALVE allows to download large quantities of product through a reciprocating movement of 180°, while the VR8V™ allows the continuous feeding of the product inside the "Y" preventing the product from being subjected to de-mixing. To avoid accumulation of the product inside the pipeline capacitive sensors can be installed for the detection of the minimum and maximum level of the product inside the pipeline. While in order to avoid that the product creates a bridge inside the pipe it is possible to install a pneumatic vibrator. On the two outlets of the "Y" you can install a DOUBLE VALVE, an SRF, a ROTARY VALVE or a VACUUM ROTARY VALVE, depending on the machine to discharge, the type of system which interfaces with the CO.RA. system and the dosing precision required. It is possible to install a flow diverter on this system in the case the working cycle requires to divert the product flow only to one of the pipelines while closing the other pipeline at the same time.*

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#### 4. DOSING SYSTEM FOR CAPSULES AND TABLETS

*The dosing system for capsules and tablets is used in processes where the the product must be dosed in bags or containers preventing the tablets from being damaged or broken during the dosing and generating waste. This type of system was conceived for all those customers that have an intermediate stocking of the product preventing it from deteriorating. The fundamental component of this system is the ROTARY VALVE TABLET and the tablet discharge system which can be customized and adapted to the customer's needs. The system can fill single bags or various bags inside of a box guaranteeing the uniformity of the weight discharged for each bag. An HMI operator panel with all the settings related to the product to discharge, counting of the number of tablets rejected and number of bags and number of containers filled. Furthermore it is possible to install a printer to print an adhesive label where, apart from all of the above indicated data, it is also possible to include the customer order number, batch number, etc... in order to maintain a continuous traceability of the product supplied to the customer as well as the quantity supplied. This system is also designed to close seal the bags and the box in a continuous manner, going from the bag sealing station to the box sealing station by means of a conveyor, which not only prevents the operator from the fatigue of having to move the containers, it also allows him to regulate the height by means of the HMI panel. This system can be equipped with air blowers for products which are particularly sensitive to humidity or a suction system which reduces the the dust deriving from the tablets to a minimum and therefore reduces the operator's contact to the product.*

Dettaglio ROTARY VALVE TABLET con tappi in silicone  
*Detail of ROTARY VALVE TABLET with silicon caps*

Dettaglio ROTARY VALVE TABLET con capsule  
*Detail of ROTARY VALVE TABLET with capsules*

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## DOSING VALVES

*Depending on the type of dosing to be executed, the following valve can be installed:*

### 1) ROTARY VALVE: For bulk dosing

*This type of valve is used when the powder or granules to discharge do not require a high level of accuracy. With this type of valve the dosing accuracy is  $\pm 500$  g.*

### 2) ROTARY VALVE TABLET:

#### *Dosing of tablets, capsules and silicone caps*

*This type of valve is used when the product to dose is in tablets or capsules and therefore subject to breaking. Thanks to its silicone rotor, this valve allows such dosing without problems.*

### 3) DOUBLE VALVE, SRF: Fine dosing

*The DOUBLE VALVE is used when the product to discharge requires a high dosing accuracy. The DOUBLE VALVE is composed of two intercepting elements, the butterfly for bulk dosing and quick discharge of a large quantity, while the rotor trickle feeds and discharges small quantities of product with an accuracy of  $\pm 50$ g.*

*The SRF is used when the product to discharge requires a high level of accuracy. The SRF is composed of two intercepting elements, the main rotor which allows quick discharge of large quantities and the smaller rotor which trickle feeds and discharges small quantities of product with an accuracy of  $\pm 50$ g*

### 4) VR8V: Continuous dosing

*The VR8V is used when the product to discharge requires a high level of accuracy, even with products that have de-mixing problems.*

*The VR8V has a spherical design which allows the accurate control of the discharge volume.*

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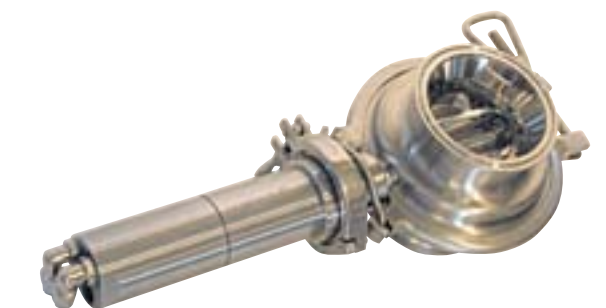
## **TYPE OF SCALE WHICH CAN BE INSTALLED**

---

Based on the type of product, on the customer's specific requests and the design of our technical dept. a scale is chosen and sized for use in a specific application.

The weighing terminal is composed of the following parts:

**Scale plate:** stainless steel structure conceived to allow cleaning so that there is no contamination or corrosion. The stainless steel platform is hermetically sealed (IP68) and is indicated for use in hazardous areas.

**Scale control panel:** the backlit LCD display quickly and clearly shows the resulting weight up to 25 mm high digits. The introduction of the data is simplified thanks to the logic of the function keys. Apart from the weight, the terminal can perform other additional important functions, such as automatic tare, data storage, functionality testing of the instrument, identification of the product and date / time function that simplify and improve operating modes.

## **TYPE OF CONTROL PANEL**

---

**Control Panel:** All the buttons necessary to the operator to dose are on the control panel of the DOSING SYSTEM along with the indicators that allow him to verify/control the condition of the machine during the working phases. The control panel of the scale which displays the discharge weight is also installed on the control panel. In the management of complex systems where the DOSING SYSTEM must be electrically interfaced with the other machines, CO.RA. uses APLC technology with Touch Screen for a simpler and more straight forward interface. The use of the Touch Screen panel offers numerous advantages compared to traditional systems such as: personalization of the system, expansion of the functionality of the machine, setting of the working cycle, batch management, use of the machine by means of a password, printing adhesive labels and all of the features that are required by the end customer.

## **DOSING SYSTEM SUPPORT**

---

There are various types of DOSING SYSTEM: mobile on wheels which allows it to be installed and/or removed when necessary for cleaning and maintaining that, depending on the work area where the operation is carried out, could cause the contamination of the environment. Furthermore, this feature allows the system to be used in various rooms. Once the system is correctly installed in its service position it is ready to be used.

## **MACHINE SAFETY**

---

The dosing system is designed to prevent accidents or dangerous situations during the use of the machine.

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## ADDITIONAL COMPONENTS OF THE DOSING SYSTEM

---

- *Machine to be served interface via derivation of the control panel of the Dosing System*
- *Air blower for the dosing of products with tendency to humidity*
- *Suction during the dosing of products with tendency to create dust*
- *Connections for inert gases*
- *Operation of system via the HMI panel*
- *Operation of system via PLC interface*
- *Sampler*

## ACCESSORIES

---

CO.RA. not only designs and produces the Dosing System, it also designs and manufactures a wide range of accessories for the system, such as :

- *Discharge pipelines*
- *Containers*
- *Loading hoppers*
- *Conveyors for the transportation of the container to the drum/container stocking area*
- *Trolleys for the transportation of drums/containers*
- *Lifting systems for bags/containers to avoid operator fatigue*
- *Bridge-breaker for products which tend to create bridges*
- *Through wall connection in case the machine to discharge is installed on the upper floor*
- *Silicone hose*
- *Sil caps*

*That adapt to the customer's specific requirements*

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---

The documentation supplied is:

- Material certificates of the parts in contact with the product
- O & M Manual
- Supply layout
- Electric and pneumatic diagrams
- P & ID
- CE certificate of conformity
- ATEX certificate of conformity

Documentation which can be supplied on request:

- Functional specifications
- FAT e SAT protocols
- IQ ed OQ protocols
- Surface finish certificate for the parts in contact with the product

The DOSING SYSTEM is in accordance with the requests of Machine Directives and can also be supplied for use in ATEX environments : zone 1/21 internally and externally

### Services co-related to the supply

- Installation
- Start-up assistance from our specialized technicians
- Personnel training
- Annual maintenance by one of our specialized technicians
- Assistance within 48 hours from your claim

The DOSING SYSTEM is designed and produced respecting the customer's requirements and to better guarantee this service it is possible to carry out tests in our facilities with the DOSING SYSTEM installed in our Technology Center.

## USEFUL INFORMATION

---

For further information please contact our Sales Dept by phone +39 0583 20590 or by e-mail: [sales@coraitaly.net](mailto:sales@coraitaly.net)

UNI EN ISO 9001:2008 - Cert. No. 5762/1

REG. No. IT-71044

ATEX 94/09/CE

CO.RA. s.r.l. - 55011 Spianate - Altopascio (Lucca) - Italy - Loc. Chiappini, 51  
Tel. +39 0583 20590 r.a. - Fax +39 0583 20481 - info@coraitaly.net - www.coraitaly.net