ROTARY
VALVE

CO,RA[®]

ROTARY VALVE

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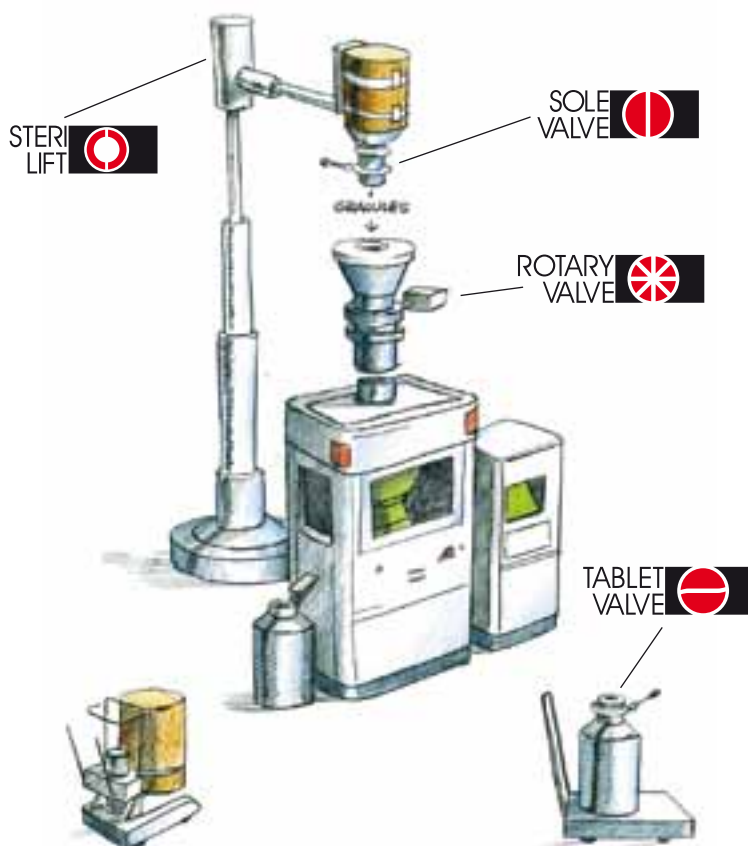
*The CO.RA.® ROTARY VALVE solves the problem of interception and control of product flows in a chemical and pharmaceutical plant.
In particular allow for machine feeding with granules and powders by equally dosing the product.*

*Its implementation derived from careful observation by CO.RA.® of the operator's requirements during the most delicate pharmaceutical processes in a sterile room, especially the cleaning phase.
The ergonomics of the parts of which it is composed make it easier for the operator during the disassembly and re-assembly of the valve.
The valve can be used in the pharmaceutical, chemical, food and cosmetics industries as a loading or discharging valve on machine or bottom valve for powder containers.*


MACCHINE SERVITE / EQUIPMENT SERVED:

- MULINI - MILLS
- COMPRIMITRICI - TABLET PRESS
- OPERCOLATRICI - FEEDING MACHINES
- MICRODOSATRICI - MICRO DOSER
- SETACCIATRICI - SIEVES

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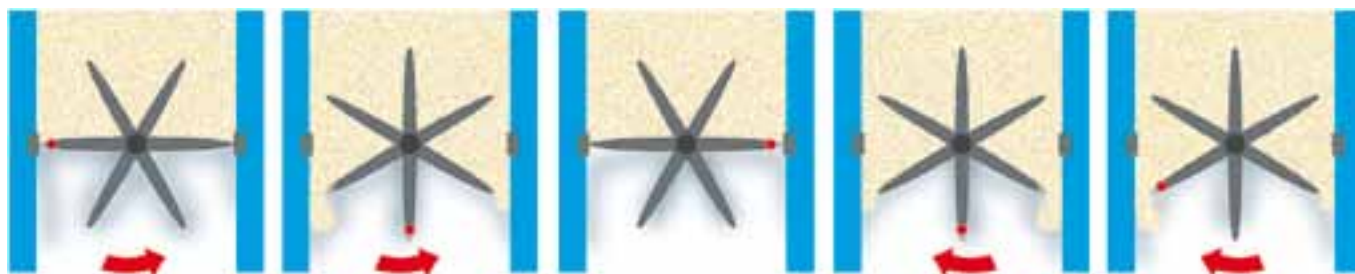Thanks to its stellar design and its movement, (alternating 180° in one direction and 180° in the other) the rotor gathers the product and transfers it to the underlying machine while simultaneously stopping the discharge of product from the upstream hopper thus avoiding the de-mixing of the product thanks to the alternate rotation.

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











Speed and stop motion can be adjusted as needed through a control panel interfaced to the valve.

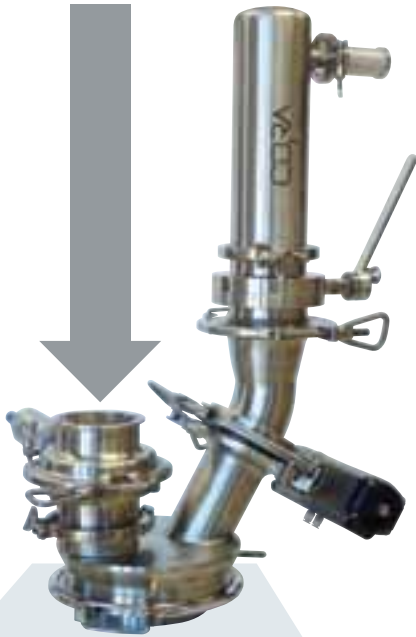The range of halfbodies available and interchangeable, allows the user to make the application more suitable according to its disassembly needs.

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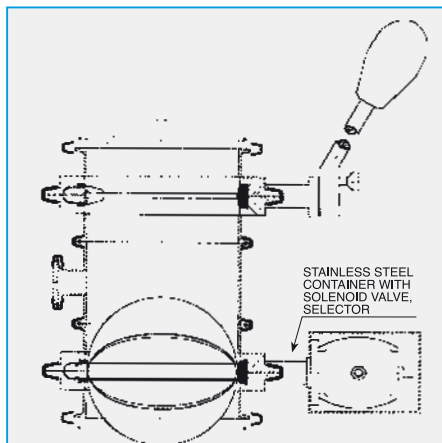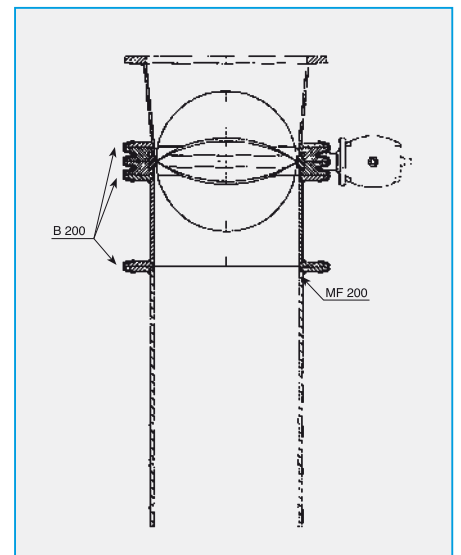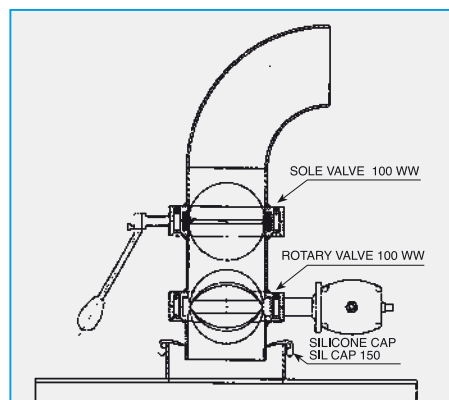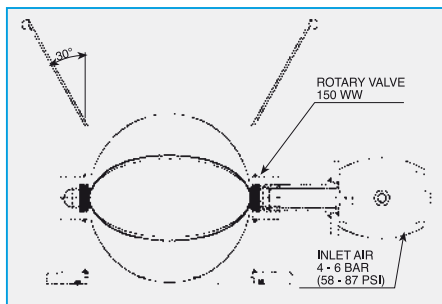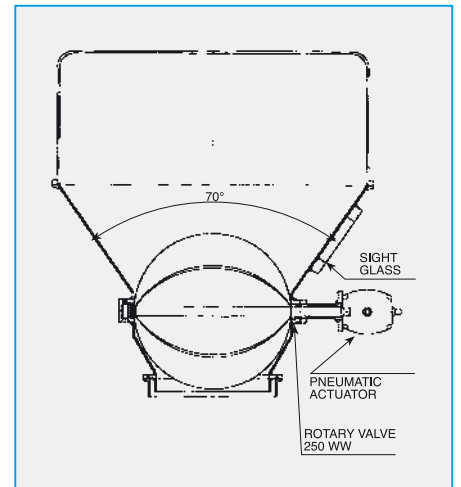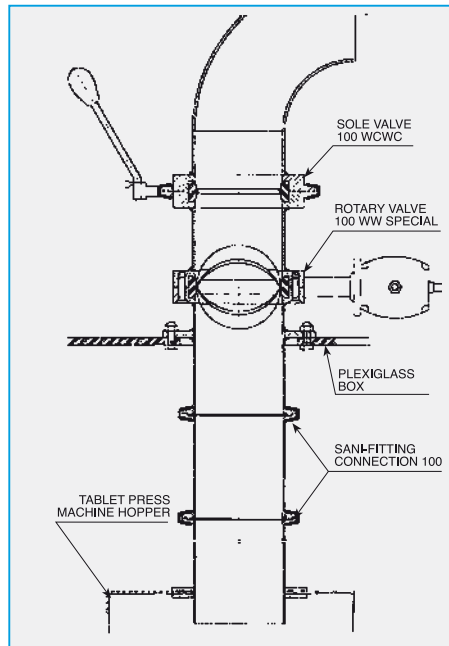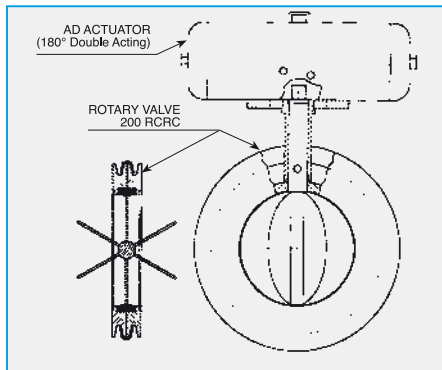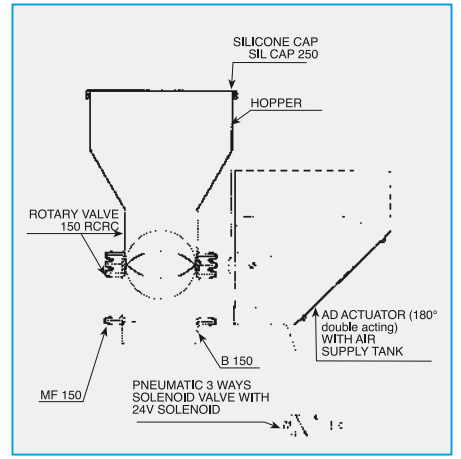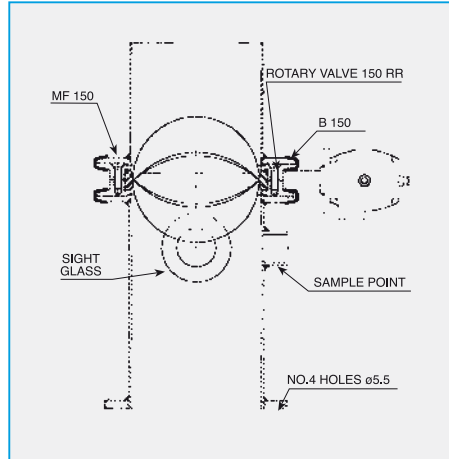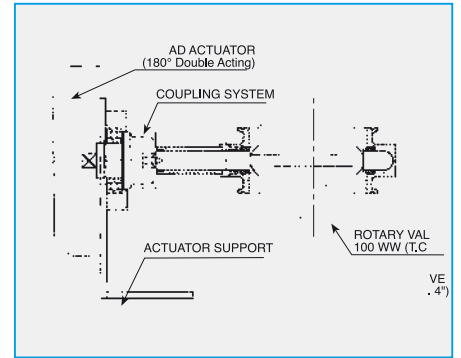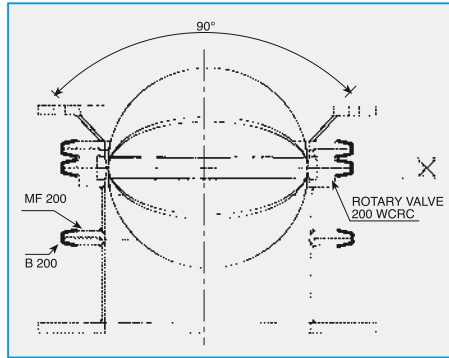- AISI 316L (DIN1.4404) satin finished body (mirror polished upon request)
- AISI 316L (DIN1.4404) mirror polished rotor
- 6 pockets (8 upon request) rotor operated by an alternating movement to feed uniform quantity of product
- No O-Ring (OR) on the shaft
- Platinum cured silicone gasket, black silicone, VITON, EPDM, PTFE core with white or black silicone, gasket with inox mesh
- Gasket for ATEX version: black silicone, PTFE core with black silicone
- Easy to disassemble, inspect, clean- particularly the "quick release" series with "CLAMP"
- Very light weight and compact
- Stainless steel and silicone
- Manufactured according to G.M.P.
- Allows continuous feeding of granules and powders without obstructing product flow
- Can handle pressure and vacuum (in closed position)
- High temperature resistant
- Interchangeability with SOLE VALVE and TABLET VALVE components
- Each valve bears the CO.RA.® trademark and identification batch number (ex.: CO.RA. I/11/298/1)

TABELLA PORTATE ROTARY VALVE - CAPACITY CHART FOR ROTARY VALVE

ND	PORTATA / CAPACITY CYCLE (Litres/cycle)					TENUTA / SEALING			COPPIA MEDIA MEAN TORQUE (n/m)
	N° Lobi Pockets number	Portata per Lobo Pockets capacity	N° Lobi Pockets number	Portata per Lobo Pockets capacity	Portata per ciclo Total per cycle	Vuoto Vacuum (mm Hg)	Pressione / Pressure		
							(bar)	(psi)	
80	6	0,040	8	0.03	0.09	650	0,2	3	20
100	6	0.09	8	0.07	0.26	650	0,2	3	23
150	6	0.29	8	0.22	0.88	650	0,2	3	30
200	6	0.7	8	0.52	2.09	650	0,2	3	35
250	6	1.36	8	1.02	4.09	650	0,2	3	60

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R = raggio del rotore(ND/2) in dm

n = numero cicli al minuto

ROTARY VALVE capacity calculation (Litres/Minute)

To calculate the capacity multiply the atmospheric volume (see plan "Total per cycle") for the cycles in a minute

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Dimensions and technical details can be subject to variation without notice
The measurements for G and L are related to D1, for dimensions of D2 add 15 mm to D1

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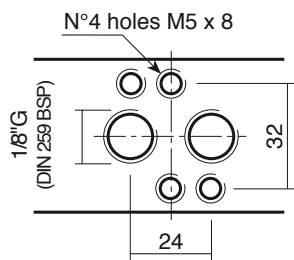
Pneumatic actuator is a driving member through which the movement is transmitted to the rotor.
The double acting actuator supplied (D.E.) drives either the opening rotation or closing rotation.
Feeding: filtrate compressed air, better if lubricated, working pressure for right working of the valves min. 5 Bar max. 6 bar.

- Upon request the ROTARY VALVE can be equipped with:
- Double Acting actuator in stainless steel carter
 - NAMUR mono or bistabile solenoid 3/2, 5/2 and 5/3 – standard or Ex rated (explosion proof)
 - Proximity, pneumatic, electric limit switches
 - Electric and/or pneumatic actuator, motor or gear unit
 - Electro-pneumatic control panel for the regulation of opening and closing speed

For more information about accessories for CO.RA.® valve actuators please consider the Sole Valve catalogue.

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*Rotary Double acting actuator (AD) of anodized aluminum to ISO 5211
 Air supply 5 ÷ 6 bar (58 ÷ 87 psi)
 Air consumption to 5.6 bar (80 psi)
 AD 63-180 = about 0.26 liters
 AD 85-180 = about 0.55 liters*

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CONTROL PANEL

The ROTARY VALVE can be supplied with pneumatic or electro-pneumatic control panel. The control panel case can be supplied in plastic material or AISI 304.

The types of control panel that can be supplied are the following:

VP PANEL

Completely pneumatic control panel, with the possibility to adjust the cycle speed (*), by means of the flow regulator installed internally.

VSP PANEL

Completely pneumatic control panel, with the possibility to adjust the pause time (**) by means of the two timers (one for the clockwise rotation and one for the counter clockwise rotation) and to adjust the cycle speed (*), by means of the flow regulator installed internally.

VSEP PANEL

Electro-pneumatic control panel, with the possibility to adjust the pause time (**) by means of the two timers (one for the clockwise rotation and one for the counter clockwise rotation) and to adjust the cycle speed (*), by means of the flow regulator installed internally. The start and the stop of the cycle take place through an external electrical signal. The timer, that allows the valve regulation is installed on the control panel door, thus allowing the regulation by the operator.

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
The mounting flange can be supplied with male end MF/M if connected to valves and female end MF/M with OR and MF/F with OR; the same for blanking plate P/M, P/F, OR are made of silicone if requested of other elastomers.
*Ask us for the list of all hand levers, supports and opening systems.
** Also other types of coupling available: Batonnet, chute, etc.

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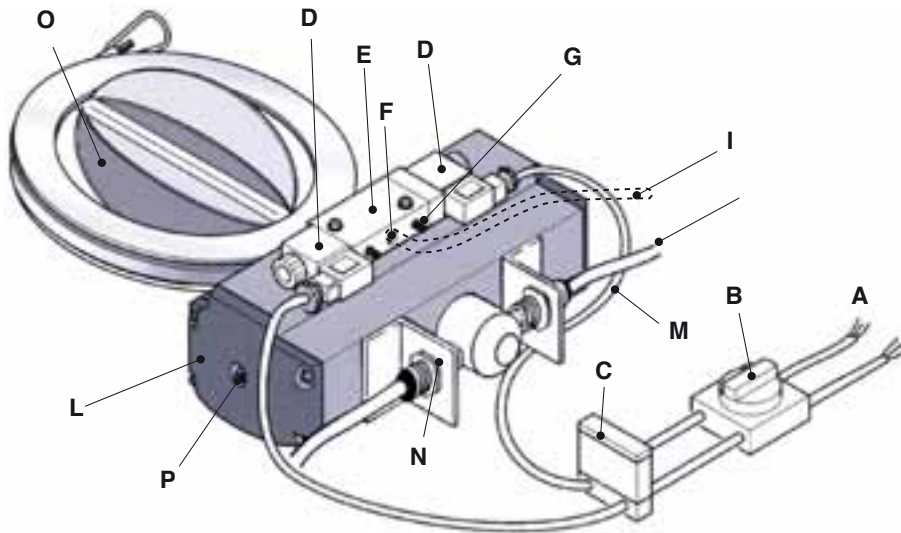
Example of assembly plan with double acting (AD) pneumatic actuator

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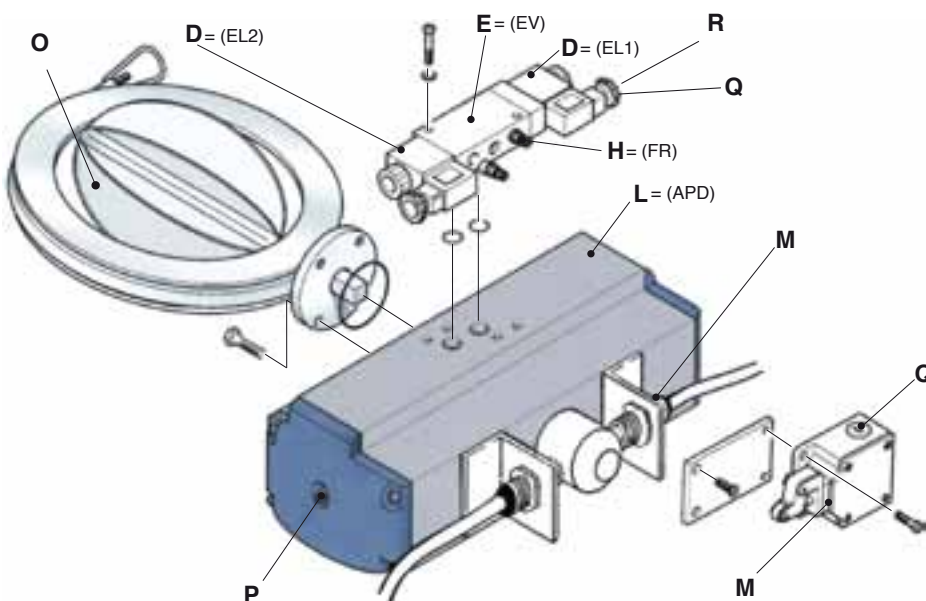
Exploded view for spare parts ordering (for spares for each main component ask for detail drawings)

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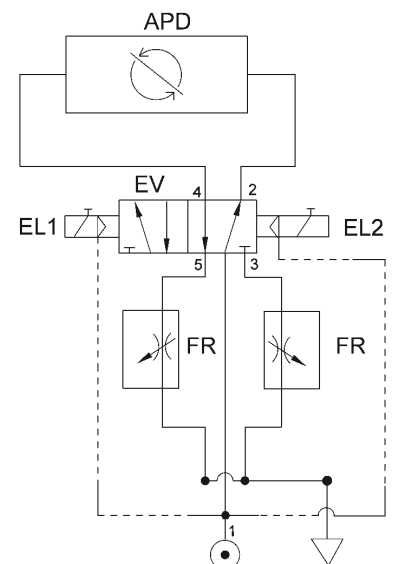DOCUMENTATION AND CERTIFICATES

The documentation that can be supplied with the valve is:

- 3.1 material certificates for the parts in contact with the product and FDA and USP certificates for the gaskets
- Operation and Maintenance Manual
- Certificate of surface roughness
- Certificate of CE conformity
- FAT and SAT protocols
- IQ and OQ protocol
- Other documentation on request

The ROTARY VALVE comply with Machinery Directive and are marked **CE**, they can also be supplied for ATEX zone. The valve markings can be as follows:

II 1GD/2GD c 85°C (T6) 0°C Ta+40°
II 1D/2GD c 85°C (T6) 0°C Ta+40°
II 2GD/2GD c IIA 85°C (T6) 0°C Ta+40°
II 2GD/2GD c IIB 85°C (T6) 0°C Ta+40°

Services related to the supply

The services related to the supply are:

- Assistance within 48 hours from receipt of mal-function report

The ROTARY VALVE are designed and manufactured respecting customer requirements and to guarantee this, it is possible to run tests in our facilities with one of the ROTARY VALVE available in our Technology Center.

USEFUL INFORMATION

We suggest you to use the above indicated abbreviations (W, WC, WR, WCRC, ...) to easy fill in your P.O.; please consider the rotor usually is made of AISI 316L polished, the supports are AISI 316L satin. The seal is elastomer.

Each valve with same diameter has interchangeability about half bodies, supports, rotors, seals, actioning, ecc.

The clamps are the same for quick clean and quick release.

Gasket available in several material and types are in accordance with FDA, for more information please ask for SOLE VALVE catalogues.

For further information please contact our Sales Dept by phone +39 0583 20590 or by e-mail: sales@coraitaly.net

Reproduction of part or all of the contents in this publication without the written consent of CO.RA. is forbidden

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*Dosing system
for plastic caps*

The ROTARY VALVE TABLET, is a special version of the ROTARY VALVE. This type of valve has been designed to solve the problem of interception, flow control or dosing of: tablets, capsules, silicone plugs, etc..

In particular it allows the continuous feeding of machines such as tablet coating machines, blister machines, intermediate filling systems, etc..

The main features of the ROTARY VALVE TABLET is the silicone rotor, which does not damage the product during discharge. There are three versions of rotor, the first two versions are characterized by a coating ("Coated B") which provides a smoother and more slippery surface than the first two versions.

The rotors are manufactured in a clean room in accordance with the existing FDA and GMP regulations.

The halfbodies as well as the weight and documentation supplied are the same as for the other valves.

ATEX CERTIFICATION

The ROTARY TABLET VALVE complies with Machinery Directive, is CE marked and also can be supplied for ATEX zones. The valve markings can be as follows:

II 1GD/2GD c 85°C (T6) 0°C Ta+40°
II 1D/2GD c 85°C (T6) 0°C Ta+40°

The ROTARY VALVE TABLET to be marked / used in ATEX environment 1GD/2GD must have the butterfly made of black conductive silicone (F.D.A. approved). For the ATEX marking 1D/2GD contact our sales department to know the available elastomers (all F.D.A. approved)

The ROTARY VALVE TABLET is designed and manufactured respecting customer requirements, and to better guarantee this service it is possible to make test at our premises with one of our ROTARY VALVE TABLET in our Technology Center.

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